

Get the most from your Kemio

Palintest

Water Analysis Technologies

Kemio

Kemio Instruction Manual



Kemio is the next generation measurement platform which can be tailored to suit your application needs.

- Enrich your test results.
- Simplify your decision making with pass/fail results.
- Protect your data.

Download the full instrument manual from the product page at www.palintest.com

Register your product to receive the latest software updates.



- Palintest



- Palintest



- PalintestMarketing

www.palintest.com

Palintest

Water Analysis Technologies

Kemio Quick Start Guide




ZIINSTKEMDIS

Register your instrument

Register your Kemio instrument online to receive the latest software, instructions and updates from Palintest. Scan the QR code or visit <https://www.palintest.com/en/product-registration>.




- 1 Fill out the required fields in the form.
- 2 Enter your instrument's serial number from the bottom of the instrument.
- 3 (Optional) check the boxes to sign up to receive the latest news from Palintest and updates about instrument servicing.
- 4 Press 'Register your product' to submit the form.

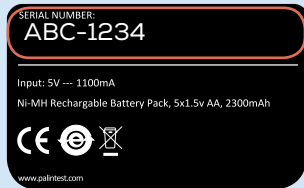

Select test types

- 1 Press CONFIGURATION, then SELECT TEST.

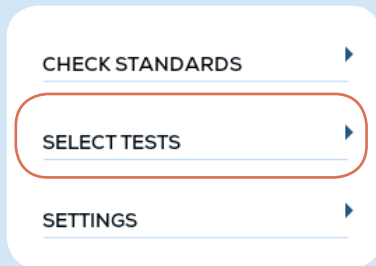

- 2 Select the test types you wish to use.



- 3 Press < to return to the CONFIGURATION menu.

Add batch information

- 1 Press BATCH INFORMATION.
- 2 Add new batch information.



- 3 Scan the QR code on the side of the sensor box.



- 4 Press  to return to the home screen.

Perform a test



Press the power button to turn on the instrument.



Press PERFORM TEST.



Follow the on-screen test instructions.



After the countdown, view the results on screen.

For more details and troubleshooting, download the full test booklet from the Palintest website at: www.palintest.com